

Scitara DLX Connector for Waters Empower 3 v2.0.0 Release Notes

Last Modified on 01/05/2024 1:54 pm EST

In this article:

- [Waters™ Empower™ 3 Connector Enhancements](#)

Waters™ Empower™ 3 Connector Enhancements

The following enhancements have been made to the Empower 3 Connector:

New Action: Retrieve Sample Set Results

The Connector now provides the ability to retrieve all results for a given sample set.

Renamed Actions

This Connector now supports the following actions:

- **Create Sample Set Method:** Creates a sample set method in Empower.
- **Retrieve Results:** Retrieves all results for the given result IDs.
- **Retrieve Result Set Results:** Retrieves all results in a result set.
- **Retrieve Sample Set Results:** Retrieves all results for a given sample set.
- **Update Sample Set Method:** Updates specific fields of an existing sample set method using either the sample name or line number as the key.

Empower 3 connector



Empower 3

This application connector is used with the Waters™ Empower™3 system to create and update sample sets and retrieve results

Version
2.0.0 (current) +

Actions (5)

Create sample set method

Creates a sample set method in Empower

Retrieve results

Retrieves results for given result IDs

Retrieve result set results

Retrieves results for given result set

Retrieve sample set results

Retrieves results for given sample set

Update sample set method

Updates specific fields of an existing sample set method using either sample name or line number as the key

Close

For more information or assistance, please contact [Scitara Service](#) to submit a ticket or use the chat bot.

© 2024 Scitara Corporation. All rights reserved. Scitara™ is the trademark of Scitara Corporation. Other trademarks mentioned are owned by their respective companies and are not associated with or sponsored by Scitara Corporation.