

DLX[™] Web Console LabX Orchestration Script v1

May 28, 2024

Module	1:	Creating	a	LabX	Orchestration
--------	----	----------	---	------	---------------

Screen	Title	On Screen	Video Time Stamp	Voice Over	SME Feedback
1.	Creating an Orchestration with the LabX Connector	Creating a LabX Orchestration		This video will demonstrate how to create a LabX Orchestration using the "Get a list of tasks" and the "Get task results" Connection actions.	
2.	Access Requirements	Access Requirements Ensure that: • A Windows File System Connection has been created • The Windows File System Connection is online		This Orchestration will include Connection actions, so ensure that a LabX Connection has been created and is online.	
3.	Creating an Orchestration with the LabX Connector: A Tutorial	Creating an Orchestration with the LabX Connector: A Tutorial		Now we will demonstrate the complete process of creating an Orchestration using the LabX Connection that we created and configured in the previous lesson. Follow the instructions and we will have a working Orchestration.	
		Video file Creating_a_LabX_Orchestration.mp4			
4.		Scitoro DXXRAre Import Import Import Import Import Import Import Import Import Import		From the DLX Orchestrations page, click Add New to create the Orchestration.	

Screen	Title	On Screen	Video Time Stamp	Voice Over	SME Feedback
5.		Scitoro LabX_Academy I Orchestration Settings Connection event ~ O Settings		Enter a name for the Orchestration.	
6.		Enabled SAVE CANCEL		Enable the Orchestration.	
7.		Scitoro		Use the trigger type, "User Triggered."	

Screen	Title	On Screen	Video Time Stamp	Voice Over	SME Feedback
8.		Scitoro		Click the plus sign and add a Connection Action step.	
9.		Connection Action Complete configuration LabX v 1.0.1(current) ~ Select a connection Select an action		Select the LabX Connector.	

Screen	Title	On Screen	Video Time Stamp	Voice Over	SME Feedback
10.		Connection Action Complete configuration LabX v1.0.1(current) ~ LabX_Academy v1.0.1 Select an action		Select the Connection we created in the previous exercise.	
11.		Connection Action Configured, check options LabX v1.0.1(current) ~ LabX_Academy v1.0.1 Get a list of tasks Returns a list of tasks created up to 30 days before the specified date		Select the "Get a list of tasks" Connection action.	
12.		Connection Action Options Options Cre Configured, check options Of 2 required options not set Cre To Date * To O5/15/2024 ···· + Status Of Task* Tring Select ···· +		In the Options tab, enter the "To Date" value in the string.	

Screen	Title	On Screen	Video Time Stamp	Voice Over	SME Feedback
13.		Connection Action Configured check options To Date* string Status Of Task* tring All XXXXXXXXXXXXXXXXXXXXXXXXXXXXXXXXXXX		Use the drop-down arrow to select the "Status of Task" value.	
14.		Enabled SAVE CANCEL		Click Save.	
15.		Signature and Reasons for change Required		Click Sign after selecting reason to complete sign off. We will then be prompted to input the username and then be redirected to the Identity Provider. (Please note that an eSign prompt will only appear for GxP implementations of DLX).	
16.		LabX_Academy Orchestration Settings User triggered ~ · · · · · · · · · · · · · · · · · ·		Click Trigger Orchestration.	

Screen	Title	On Screen	Video Time Stamp	Voice Over	SME Feedback
17.		 LabX Academy # Started by Sobiel Ansair less than a minute ago, completed in less than 30 seconds Orchestration execution completed I User Triggered ther-friggered ther-friggered Control Action Certifies action compared Cold Bings table T1967 501-502-502-5025-0055000b¹ indemnal T1967 control T1967 501-5025-0055000b¹ indemnal T1967 control T1967 504-5025-0055000b¹ indemnal T1967 control T1967 504-5025-0055-0055000b¹ indemnal T1967 control T1967	6	Expand the Orchestration execution, then select the "Get a list of tasks" Connection Action and click the drop-down arrows to view the step output data and verify execution.	
18.		scitture Sate Image: Scitture Sate	And Ansani	Click Close .	

Screen	Title	On Screen	Video Time Stamp	Voice Over	SME Feedback
19.		Scitoro		Click the plus sign to add another Connection Action step.	
20.		Connection Action Complete configuration LabX v10.1(current) ~ Select a connection		Select the LabX Connector.	

7 | Page

Screen	Title	On Screen	Video Time Stamp	Voice Over	SME Feedback
21.		Connection Action Complete configuration LabX v10.1(current) ~ LabX_Academy v10.1 Select an action		Select the Connection we created in the previous exercise.	
22.		Connection Action Configured, check options LabX v1.0.1(current) ~ LabX_Academy v1.0.1 Get task results Get the results for a task		Select the "Get task results" Connection action.	

Screen	Title	On Screen	Video Time Stamp	Voice Over	SME Feedback
23.		Connection Action Configured, check options Task internal ID thring Query Tasks From thring To Enter the text + Task GUID thring To Enter the text + Task GUID thring To Enter the text + To Enter the text +		In the Options tab, change the Task GUID expression data type to a Reference.	

Screen	Title	On Screen	Video Time	Voice Over	SME Feedback
			stamp		
24.		Select a reference value Image: Construction of the second s		To select a reference value, expand "Steps" and drill down to "result 1." Click the download arrow icon next to "taskID".	
25.		Enabled SAVE CANCEL		Click Save.	
26.		Signature and Reasons for change Required		Click Sign after selecting reason to complete sign off. We will then be prompted to input the username and then be redirected to the Identity Provider. (Please note that an eSign prompt will only appear for GxP implementations of DLX).	

Screen	Title	On Screen	Video Time Stamp	Voice Over	SME Feedback
27.		User triggered Trigger orchestration Cancel Trigger Crchestration Cancel Trigger Crchestration		Trigger the Orchestration.	
28.		Schermannersensensensensensensensensensensensensens		Expand the Orchestration execution, then select the "Get Task Results" Connection Action and click the drop-down arrows to view the step output data and verify execution.	

Screen	Title	On Screen	Video Time Stamp	Voice Over	SME Feedback
29.		Scherard		Click Close .	
30.		Sciture Sciture Sciture Science Scienc		Close the Orchestration then view the Orchestrations Execution window to verify execution.	
31.		Add Filter O UI Jäkos results () Jäkos o O 13 Jäkos o O 10		Click on the Orchestration execution then expand the "Get Task Results" Connection Action. Click the drop-down arrows under "samples" to view the name for "sampleID1", then expand "scalar" and view the data.	

12 | Page

Screen	Title	On Screen		Video Time	Voice Over	SME Feedback
				Stamp		
32.		Global Parameters	● — □ ×		Open the LabX application and double click the completed task. Verify that the Sample ID matches the data in the Orchestration execution.	
		Marked In Linne				
		metro in more				
		Method name Virtual Weight				
		Task name SW Test2				
		Tesk comment				
		Treat as priority task				
		Centinopus run				
		Scope 1 Sample scope				
		Number of samples 1				
		+ Ant @ Bernes + Mounths - L Moun Down - W length	C - M Zuccer			
		No. Sample ID				
		5W0002				
			Canopl			
33.	Results Q. Search text. Drag a column fr Creation time 26-04-2024 18: 26-04-2024 18: 26-04-2024 18: 26-04-2024 18:	Results			that the Sample ID for the task	
		Ne service tett.				
		Drag a column header here to group by that column			matches the data in the	
		Creation time Task Sam	nple ID 1 Sample		Orchestration execution.	
		26-04-2024 18:18:32 SW Test3 SW	0003 1			
		26-04-2024 18:12:19 SW Test2	1 1000			
		26-04-2024 18:06:15 T192_Flow Rate 123	1			
		26-04-2024 18:02:54 SW Test1 SW0	0001 1			
		07.11.2022 11.07/0 T171 Simela Malekine (with Result) S.	4			

Screen	Title	On Screen	Video Time Stamp	Voice Over	SME Feedback
34.		X Andre dogl, Handradan - 0 X Hand Andre dogl, Handradan - 0 X Handrad Andre dogl, Handradan - 0 X Handrad - 0 X - 0 X Handrad - 0 X - 0 X - 0 X Handrad - 0 X		Double click the task to view the details. Verify that the "Results Net Weight" value matches the data in the Orchestration Execution.	
35.		A Recap: Creating an Orchestration with the LabX Connector		We have successfully created a LabX Orchestration using the "Get a list of tasks" and the "Get task results" Connection actions.	
36.		In case of any queries,⊥please contact us at support@scitara.io		Thanks for watching.	