




Equipment: Air Clean 600 PCR Work Station

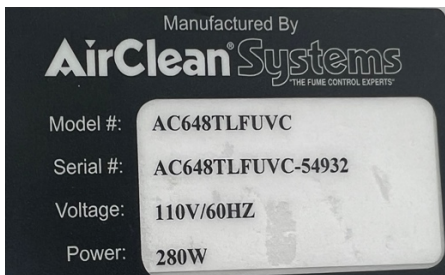

Service Line: 800-849-0472, service@aircleansystems.com

Maintenance:

1. Monthly pre-filter check for dust accumulation. 2 pre-filters located on top of the hood. Change filters every 3 months or when visible dust begins to accumulate.



TOP OF FILTER



BOTTOM OF FILTER



2 PRE-FILTERS ON TOP OF PCR HOOD



2. Change HEPA filter every 2 years. First filter was put in place 12.05.21. Next change is due by 12.04.24.
3. HEPA filter code input. As of 05.18.23 a glitch occurred with the HEPA filter code. Manually entered a 1 year code provided by them. Filter timer valid till 05.16.24. Alarm will go off once code expires. Call the service line 800-849-0472 and explain this situation and request a code for 6 months when this happens. The info. on our equipment should have a note attached on their end regarding this.
4. During a power outage, power surge, machine rebooting, etc. HEPA code may get erased resulting in a “false alarm” indicating an expired HEPA filter. Call service line and explain situation and get a new code.
5. UV Lamp Alarm will go off after 900 hours of usage to signal replacement needed. (~Early 2025)
6. Check time remaining for HEPA filter and UV lamp by checking the screen on the PCR hood. The mode button allows you to switch between the different settings. **Check timer and order replacements BEFORE time runs out and alarm starts blaring!!!**

